



pTagCFP-C vector

The vector sequence has been compiled using the information from sequence databases, published literature, and other sources, together with partial sequences obtained by Evrogen. This vector has not been completely sequenced.

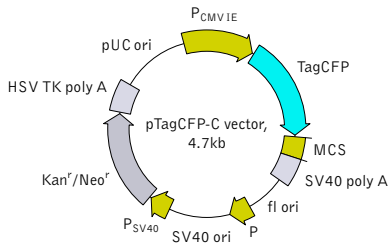

For vector sequence, please visit our Web site at <http://www.evrogen.com/products/vectors.shtml>

Multiple cloning site (MCS)

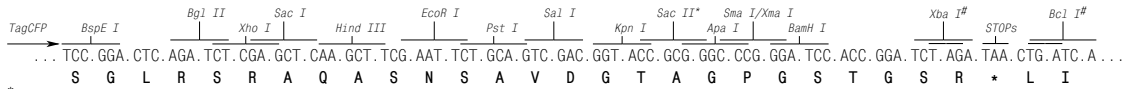

* — not unique sites.

— sites are blocked by *dam* methylation. If you wish to digest the vector with these enzymes, you will need to transform the vector into a *dam*⁻ host and make fresh DNA.

Location of features

P_{CMV IE}: 1-589
 Enhancer region: 59-465
 TATA box: 554-560
 Transcription start point: 583
 TagCFP
 Kozak consensus translation initiation site: 606-616
 Start codon (ATG): 613-615; Stop codon: 1405-1407
 Last amino acid in TagCFP: 1324-1326
 MCS: 1327-1404
 SV40 early mRNA polyadenylation signal
 Polyadenylation signals: 1547-1552 & 1576-1581
 mRNA 3' ends: 1585 & 1597
 f1 single-strand DNA origin: 1644-2099
 Eukaryotic promoter for expression of Kan^r gene
 -35 region: 2161-2166; -10 region: 2184-2189
 Transcription start point: 2196
 SV40 origin of replication: 2440-2575
 SV40 early promoter
 Enhancer (72-bp tandem repeats): 2273-2344 & 2345-2416
 21-bp repeats: 2420-2440, 2441-2461 & 2463-2483
 Early promoter element: 2496-2502
 Major transcription start points: 2492, 2530, 2536 & 2541
 Kanamycin/neomycin resistance gene
 Neomycin phosphotransferase coding sequences:
 Start codon (ATG): 2624-2626; Stop codon: 3416-3418
 G->A mutation to remove Pst I site: 2806
 C->A (Arg to Ser) mutation to remove BssH II site: 3152
 Herpes simplex virus (HSV) thymidine kinase (TK) polyadenylation signal
 Polyadenylation signals: 3654-3659 & 3667-3672
 pUC plasmid replication origin: 4003-4646

References

Gorman, C. (1985). "High efficiency gene transfer into mammalian cells." In: *DNA cloning: A Practical Approach, Vol. II*. Ed. by Glover. (IRL Press, Oxford, U.K.) Pp. 143–190.

Haas, J. et al. (1996) "Codon usage limitation in the expression of HIV-1 envelope glycoprotein." *Curr Biol*, 6 (3): 315–324 / pmid: 8805248

Kozak, M. (1987) "An analysis of 5'-noncoding sequences from 699 vertebrate messenger RNAs." *Nucleic Acids Res*, 15 (20): 8125–8148 / pmid: 3313277

Product	Cat.#	Size
pTagCFP-C vector	FP111	20 µg
Vector type	mammalian expression vector	
Reporter	TagCFP	
Reporter codon usage	mammalian	
Promoter for TagCFP	P _{CMV IE}	
Host cells	mammalian	
Selection	prokaryotic - kanamycin eukaryotic - neomycin (G418)	
Replication	prokaryotic - pUC ori eukaryotic - SV40 ori	
Use	TagCFP expression in mammalian cells; generation of fusions to the TagCFP C-terminus	

Vector description

pTagCFP-C is a mammalian expression vector encoding cyan fluorescent protein TagCFP. The vector allows generation of fusions to the TagCFP C-terminus and expression of TagCFP fusions or TagCFP alone in eukaryotic (mammalian) cells.

TagCFP codon usage is optimized for high expression in mammalian cells (humanized) [Haas et al. 1996]. To increase mRNA translation efficiency, Kozak consensus translation initiation site is generated upstream of the TagCFP coding sequence [Kozak 1987]. Multiple cloning site (MCS) is located between TagCFP coding sequence and SV40 polyadenylation signal (SV40 polyA).

The vector backbone contains immediate early promoter of cytomegalovirus (P_{CMV IE}) for protein expression, SV40 origin for replication in mammalian cells expressing SV40 T-antigen, pUC origin of replication for propagation in *E. coli*, and f1 origin for single-stranded DNA production. SV40 polyadenylation signals (SV40 poly A) direct proper processing of the 3'-end of the reporter mRNA.

SV40 early promoter (P_{SV40}) provides neomycin resistance gene (Neo^r) expression to select stably transfected eukaryotic cells using G418. Bacterial promoter (P) provides kanamycin resistance gene expression (Kan^r) in *E. coli*. Kan^r/Neo^r gene is linked with herpes simplex virus (HSV) thymidine kinase (TK) polyadenylation signals.

Generation of TagCFP fusion proteins

A localization signal or a gene of interest can be cloned into MCS of the vector. It will be expressed as a fusion to the TagCFP C-terminus when inserted in the same reading frame as TagCFP and no in-frame stop codons are present. TagCFP-tagged fusions retain fluorescent properties of the native protein allowing fusion localization *in vivo*. Unmodified vector will express TagCFP when transfected into eukaryotic (mammalian) cells.

Note: The plasmid DNA was isolated from *dam*⁺-methylated *E. coli*. Therefore some restriction sites are blocked by methylation. If you wish to digest the vector using such sites you will need to transform the vector into a *dam*⁻ host and make fresh DNA.

Expression in mammalian cells

pTagCFP-C vector can be transfected into mammalian cells by any known transfection method. CMV promoter provides strong, constitutive expression of TagCFP or its fusions in eukaryotic cells. If required, stable transformants can be selected using G418 [Gorman 1985].

Propagation in *E. coli*

Suitable host strains for propagation in *E. coli* include DH5alpha, HB101, XL1-Blue, and other general purpose strains. Plasmid incompatibility group is pMB1/ColE1. The vector confers resistance to kanamycin (30 µg/ml) to *E. coli* hosts. Copy number in *E. coli* is about 500.

Notice to Purchaser:

TagCFP-related materials (also referred to as "Products") are intended for research use only. The Products are covered by European Pat. 06809023 and other Evrogen Patents and/or Patent applications pending. By use of these Products, you accept the terms and conditions of the applicable Limited Use Label License #001: <http://www.evrogen.com/products/Evrogen-FP-license.shtml>. The CMV promoter is covered under U.S. Patents 5,168,062 and 5,385,839, and its use is permitted for research purposes only. Any other use of the CMV promoter requires a license from the University of Iowa Research Foundation, 214 Technology Innovation Center, Iowa City, IA 52242.

MSDS information is available at <http://www.evrogen.com/MSDS.shtml>